MACHEREY-NAGEL

Crimping tools




For tight and safe sealing

- Manual crimping tools
- Electronic crimping tools
- Related products
- Troubleshooting



Please choose the most suitable crimping tool for your individual requirement profile

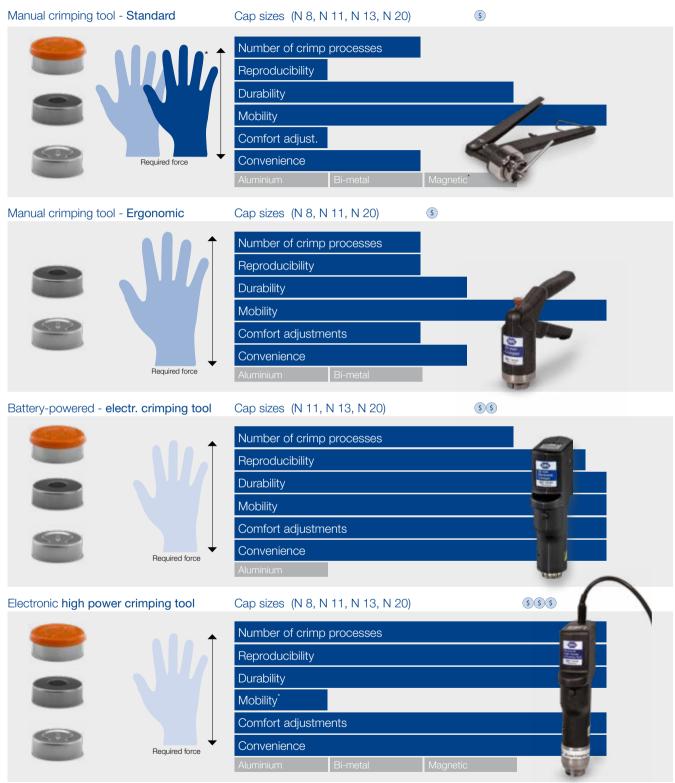

^{*} normally stationary usage with a stand

Optimal crimping tool - overview

Manual crimping tools

	Manual, ergonomic crimping tools	Manual crimping tools, standard version		
	THE STATE OF THE S	BH CONTRACTOR OF THE PARTY OF T		
Suitability / availability	For all cap materials	For all cap materials		
	For N 11 and N 20; for N 8 crimper only	For N 8, N 11, N 13 and N 20		
		For 13 mm and 20 mm Flip Top/Flip Off caps		
Sample quantities	Small series	Small series		
Criterion	One crimper and decapper each per cap size	One crimper and decapper each per cap size		
Adjustment of the crimper	Adjustment by a knob on the crimping head (easily	Crimping height adjustable in the crimping head by a hexagon key		
	accessible and well visible)	Crimping pressure adjustable by a screw on the bottom handle		
For further information see	Page 4	Page 5		
Additional information	www.mn-net.com/manualcrimper	www.mn-net.com/manualcrimper		

Electronic crimping tools

	Electronic crimping tools (battery-powered)	Electronic high power crimping tool (with fixed power supply)
Suitability / availability	Only for aluminum crimp caps N11, N13 and N20 (not suitable for magnetic/bi-metal crimp caps)	For all cap materials For N 11, N 13 and N 20; for N 8 crimping head only
	For 13 mm and 20 mm Flip Top/Flip Off caps	For 13 mm and 20 mm Flip Top/Flip Off caps
Sample quantities	Small and medium size series	Medium and large size series
Criterion	One crimping and decapping tool each per cap size	Exchangeable crimping / decapping heads
Adjustment of the crimper	Simple adjustment by pushing the up and down arrows on top of the tool (underneath the OLED display)	Simple adjustment by pushing the up and down arrows on top of the tool (underneath the OLED display)
	Settings mode with language selection, statistics, log data and different reset options	Settings mode with language selection, change of jaw set, statistics, log data and different reset options
For further information see	Page 6	Page 7
Additional information	www.mn-net.com/electroniccrimper	www.mn-net.com/electroniccrimper

For technical questions about our crimping tools and / or vials and caps you are always welcome to contact us under vials@mn-net.com.

Manual crimping tools

Ergonomic version

Features

- More lightweighted than complete steel crimpers
- Ergonomically designed handles
- Adjustment of the crimping pressure by a knob on the crimping head that is easily accessible and well visible. Crimp setting doesn't need any further fixation.
- Activated by bottom handle motion only which allows a steadier and safer hold of the tool during crimping
- Due to design and alignment of the crimping head better vertical clearance over the vial
- Brand and product identification on the front label of the tool



The adjustment knob, which sets the stroke of the tool and thus the tightness of the crimp, is conveniently positioned on top of the crimping head. The easily-viewed knob clearly displays + and - symbols along with directional arrows, in order to simplify adjusting the level of crimp desired. During the crimping process, in which the bottom handle is pulled up, the adjustment knob moves towards the crimper body. Once the knob touches the head, the crimp setting has been reached and the crimping process stops.





Additional information on adjustment and handling of the tool

Instruction leaflet as pdf download on the respective product detail section under the "Downloads" tab.

Ordering information

Pack of	REF
1	735208
1	735211
1	735220
1	735311
1	735320
1	735509
	Pack of 1 1 1 1 1 1 1

Our crimp neck vials and closures N11, N13 respectively N20 you'll find ...

...on our website under www.mn-net.com / vials



A selection of the most common products you will find on page 8.

Manual crimping tools

Standard version

Features

- Adjustable crimping height by a hexagon key, which allows to move the inner part of the crimping head up and down (not possible for manual crimpers N 8)
- Crimping pressure adjustable by a screw in the handle; fixation of the crimp setting by a nut
- Manual crimpers for N 13 and N 20 Flip Top / Flip Off caps also available
- Robust metal crimper with a long life time
- Blocky handles allow grabbing and pressing of the tool handles at any position



Correction of inaccurate crimp features

Crimp too loose (undercrimped)

→ Adjust crimping height (in the crimping head, see ②) and/or crimping pressure (by screw in the handle, see ①)

Septa emerges through the center hole (overcrimped)

→ Adjust crimping height (in the crimping head, see ②) and/or crimping pressure (by screw in the handle, see ①)

→ Adjust crimping height (in the crimping head, see ②)

→ Adjust crimping height (in the crimping head, see ②)

→ Adjust crimping pressure (by screw in the handle, see ①)

Adjustment of crimping pressure

- Further unscrew the screw ① in the handle → handles cannot be pressed so close together → lower crimping pressure
- 2. Further screw the screw ① in the handle → handles can be pressed closer together → higher crimping pressure

Adjustment of crimping height

- 1. Fixate the handles of the crimper with the striker (3).
- 2. Hold crimping head ② firmly in one hand and insert the supplied hexagon key in the screw inside the head.
- 3. a) Make a half up to a complete turn to the left (counter clockwise rotation), in case the crimp is too loose.
 - b) Turn the hexagon key to the right (clockwise rotation), in case the crimp is too tight.

Additional information on adjustment and handling of the tool

Instruction leaflet as pdf download on the respective product detail section under the "Downloads" tab.

Ordering information

Description	Pack of	REF
Manual crimpers (standard)		
Crimper for 8 mm crimp caps	1	735126
Crimper, height adjustable, for 11 mm crimp caps	1	735111
Crimper, height adjustable, for 13 mm crimp caps	1	735113
Crimper, height adjustable, for 13 mm Flip Top/Flip Off caps	1	735133
Crimper, height adjustable, for 20 mm crimp caps	1	735120
Crimper, height adjustable, for 20 mm Flip Top/Flip Off caps	1	735132
Manual decappers (standard)		
Decapper for 8 mm crimp caps	1	735408
Decapper for 11 mm crimp caps	1	735911
Decapper for 13 mm crimp caps	1	735913
Decapper for 20 mm crimp caps	1	735920

Electronic crimping tools

Electronic crimping tools (battery-powered)

Features

- Battery-powered, electronic crimping tools for 11 mm, 13 mm and 20 mm aluminum crimp caps (not suitable for 20 mm magnetic/bi-metal crimp caps; for these please see electronic high power crimping tool REF 735700 with exchangeable crimping heads)
- Mobile tools for consistent and reproducible crimp results
- OLED Display for convenient handling
- Settings mode with language selection, statistics, log data and different reset options
- Crimping pressure adjustable by pushing the up and down arrows on top of the tool
- Brushless motor for a long lifetime
- Long lasting lithium ion cell batteries (full battery charge for several hundred vials)
- Life time of battery > 1500 charges, charging time: approx. 1–2 hours; tool can be used during charging
- No serviceable parts; replacement battery separately available
- Various plug clips for charging in almost all countries come along with the tool
- CE certificate of conformity along with one year warranty
- One tool each necessary for crimping and for decapping

Adjustment and handling

Check the crimped vial for satisfactory form and tightness. If the cap spins easily, press the up arrow two or three times. Try the new setting with a new vial and cap.

Additional information on adjustment and handling of the tool

Instruction leaflet as pdf download on the respective product detail section under the "Downloads" tab.

Too Tight Too Loose

Ordering information

Description	Pack of	REF
Electronic crimpers (battery-powered)		
Electronic crimper for 11 mm aluminum crimp caps	1	735511
Electronic crimper for 13 mm aluminum crimp caps	1	735513
Electronic crimper for 13 mm Flip Top / Flip Off caps	1	735533
Electronic crimper for 20 mm aluminum crimp caps	1	735520
(not suitable for magnetic/bi-metal crimp caps)		
Electronic crimper for 20 mm Flip Top / Flip Off caps	1	735532
Electronic decappers (battery-powered)		
Electronic decapper for 11 mm aluminum crimp caps	1	735611
Electronic decapper for 13 mm aluminum crimp caps	1	735613
Electronic decapper for 20 mm aluminum crimp caps	1	735620
(not suitable for magnetic / bi-metal crimp caps)		
Accessories for battery-powered, electronic crimping and decapping tools		
Replacement battery 6.6 Volt, 8.6 Wh	1	735500
Stand for electronic crimping tools	1	735501
Rack for two electronic crimping tools (battery-powered)	1	735509



Electronic crimping tools

Electronic high power crimping tool with exchangeable crimping / decapping heads for all 11 mm, 13 mm and 20 mm crimp caps (aluminum, magnetic, bimetal)

Features

- Due to a more powerful motor suitable for crimp caps made of different materials
- With fixed power supply
- Exchangeable 11 mm, 13 mm and 20 mm crimping and decapping heads with quick-release jaw sets (need to be ordered separately)
- OLED display with language selection, change of jaw set, statistics, log data and different reset options
- Brushless motor for a long lifetime
- No serviceable parts
- CE certificate of conformity and one year warranty
- For more convenient handling a stand is optionally available



Instruction leaflet as pdf download on the respective product detail section under the "Downloads" tab.



Ordering information

0		
Description	Pack of	REF
Electronic high power crimping tool		
Electronic high power crimping tool with fixed power supply	1	735700
Accessories for electronic high power crimping tool 735700		
Crimping head for 11 mm crimp caps	1	735711
Crimping head for 13 mm crimp caps	1	735713
Crimping head for 13 mm Flip Top / Flip Off caps	1	735733
Crimping head for 20 mm crimp caps	1	735720
Crimping head for 20 mm Flip Top / Flip Off caps	1	735732
Decapping head for 11 mm crimp caps	1	735811
Decapping head for 13 mm crimp caps	1	735813
Decapping head for 20 mm crimp caps	1	735820
Stand for electronic crimping tools	1	735501

In case you are interested to see how to handle our electronic crimping tools (battery-powerded), you can watch a demo video under www.mn-net.com/electroniccrimper.

Crimp neck vials and closures · Selection ■N 11 and ■ N 20

Crimp neck vials	Is N11, wide openin	g (image scale 1:2	11 8 × 32 mm	11 & x 32 mm	1	₩ 1 (10 5 50 mm
70201HP 70201HP.2	702885 / 7020 702892 / 7020		702141	702015 702016	70:	2134	702891 702014
1.5 mL	1.5 mL	1.1 mL	1.1 mL	1.1 mL	0.2	2 mL	0.2 mL
	label + scale ¹ silanized	15 µL funn glass botto		conical, r pedestal	glass plate inte	nsp., with egrated 0.2 mL ass insert	with integrated 0.2 mL insert
Ready assemble	ed crimp closures N	l 11, aluminum, sil	ver, center hole (imag	ge scale 1:2)			
70284	702730	70256	70231	702001	70288	702823	702995
PTFE virginal, white, 0.25 mm	Red Rubber/ FEP colorless, 1.0 mm	Natural Rubber/ Butyl red-orange TEF colorless; 1.0 mm	Natural Rubber/ / Butyl red-orange/ TEF colorless; 1.3 mm	red-orange/	Silicone white/ PTFE red, 1.3 mm	Silicone white/ PTFE blue, cross-slit, 1.5 mm	PTFE red / Silicone white / PTFE red, 1.0 mm
Crimp neck vial	s N 20 (image scale	1:2)					
70206.36/7021	(_{general park}) 1.36 70254	702	Garden Ga	25.6 x 75.6 mm (DANI, Agient")	22 6 x 75 6 mm (CTC, Martar?)	02263.2	0208.36
20 mL	20 mL	20	mL	20 mL	20 mL	5	0 mL
flat bottom	rounded bo		nded bottom	flat bottom	rounded b		at bottom
flat DIN crimp ne		·	eled HS crimp neck	flat DIN crimp neck	flat DIN cri	mp neck fl	at DIN crimp neck
Ready assemble	ed crimp closures, p	olain crimp caps a	nd single septa N 20	(image scale 1:2)			
Butyl re PTFE g		utyl light gray / IFE dark gray	Molded septa Butyl	/ Butyl dark gra PTFE gray		e blue transpa- TFE colorless	Silicone white/ PTFE beige
7027	7	702057	702101	702D20TB	-	702780	70278
A	Assembled in N 20 alu	ıminum crimp caps	center hole (REF 702	804 without liner)			
70277	73	702775	70234.9	70234.10/702	234	702093	702094
A	Assembled in N 20 alu	ıminum pressure re	ease caps, center hole	e (REF 702799 withou	it liner)		
70283	36	702829	70234.8	702071	-	702927	702835
A	Assembled in N 20 bi-	metal crimp caps, o	center hole (REF 7028)	33 without liner)			
_	Accombined in MOC	702838	- contar hale (DEE 700)		7	702834	702837
A	Assembled in IV 20 M8	agnetic crimp caps,	center hole (REF 7028	ooo wiinou(iiner)			
70277	74	702928	_	702928.9	-	702929	_

All about crimping and decapping of crimp neck vials

Crimpers from MACHEREY-NAGEL are supplied with a basic adjustment for the individual cap size and the septa thickness that is typically used for these caps. Crimp necks and septa thicknesses, however, can differ from manufacturer to manufacturer. Furthermore special consideration with regard to 20 mm crimp caps needs to be paid to the fact that these are used in combination with vials having a flat DIN crimp neck as well as with those having a beveled top Headspace crimp neck (with a very low bearing surface for the septa). The different crimp neck designs require different crimp settings of the tool.

Due to the above further adjustments of the tools might be necessary, in order to achieve an optimal crimp result.

An optimal crimp result is characterized by the following features:

- 1. The cap surface is flat and the cap sides fit tightly and firmly around the glass edges without any deformation of the aluminum surface.
- 2. The septa neither emerges out of the center hole of the cap nor is being sucked into the vial.
- 3. The cap cannot be turned by usage of single fingers without enormous expenditure of energy.

Beveled Headspace crimp neck



Flat DIN crimp neck



Good crimping result





Crimp too tight







Crimp too loose





Often the so-called "twist test" of the crimp cap is being used, in order to verify the crimp result. This test is only of limited relevance. On one hand the slippery PTFE lamination of the septa results already by a relatively low expenditure of energy in a torque that enables the turn of the cap, even if the cap was crimped correctly. As the test often is done with the complete hand (instead of using single fingers) and with a high expenditure of energy, the false statement is even strengthened.

On the other hand overcrimped closures that cannot be turned, which have, however, either deformed cap sides or sucked in septa, may not be as tight as you would expect from an optimal crimp. This is due to the fact that septa material is sucked into the vial opening by the far too high crimping pressure and is then, however, missing at the cap sides for sealing. The same is also valid for septa that - due to a wrong adjustment of the crimping height and a too strong crimping pressure - emerge through the cap's center hole.

How appropriate adjustments of the crimping tools can be carried out, depends on the type of crimping tool and can be taken from the corresponding instruction leaflet.

General instructions:

- For beveled top Headspace crimp necks the crimper needs to be adjusted tighter than for flat DIN crimp necks
- Decappers don't need any adjustment

For more information on optimal crimping please ask for our poster with MN REF KATDE/EN200153.

www.mn-net.com



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6-8 · 52355 Düren · Deutschland DE/International: CH: FR:

+49 24 21 969-0 Tel.: +49 24 21 969-199 E-Mail: info@mn-net.com

+41 62 388 55 00 Tel.: Fax: +41 62 388 55 05 E-Mail: sales-ch@mn-net.com

+33 388 68 22 68 Tel.: +33 388 51 76 88 Fax: E-Mail: sales-fr@mn-net.com

Tel.: +1 484 821 1272 Fax:

